Micrococcal Nuclease

Code No. 2910A **Size:** **15,000 U**
Conc.: **20 U/ μ l**

Supplied Reagents:
10X Micrococcal Nuclease Buffer 1 ml

Description :

Micrococcal Nuclease is an endonuclease that preferentially digests single-stranded DNA and RNA especially at AT or AU rich region, but will also digests double-stranded DNA. This enzyme digests 5'-phosphodiester bonds of DNA and RNA, and yield 3'-phosphate mononucleotides and oligonucleotides. This enzyme requires Ca²⁺ and is completely inactivated by EDTA or EGTA.

Storage Buffer :

50 mM Tris-HCl, pH 8.0
1 mM DTT
50 mM NaCl
50 % Glycerol

Storage: -20°C

Source: *Staphylococcus aureus*

Unit definition :

One unit is the amount of enzyme that produce 1 OD₂₆₀ of acid-soluble products in 30 minutes at 37°C and pH 8.0, with heat-denatured calf thymus DNA as the substrate.

Reaction mixture for unit definition :

20 mM Tris-HCl, pH 8.0
5 mM NaCl
2.5 mM CaCl₂
0.5 mg/ml substrate DNA

Quality Control Data :

Please see the Certificate of Analysis (CoA) for each lot. You can download the CoA on Takara Bio website.

Applications :

1. Digestion of nucleotides included in crude cell extract.
2. Digestion of chromatin for nucleosome preparation.

Note :

This enzyme requires Ca²⁺ and completely inactivated by EDTA or EGTA.

Composition of Supplied Reagents (Store at – 20°C) :

10X Micrococcal Nuclease Buffer
200 mM Tris-HCl, pH 8.0
50 mM NaCl
25 mM CaCl₂

* The supplied buffer gives the same composition as that used for the unit definition.

References :

- 1) Heins J N, *et al. Proc Nucl Acids Res.* (1966) **1**: 79.
- 2) Pelham H R and Jackson R J. *Eur J Biochem.* (1976) **67**: 247.
- 3) Telford D J and Stewart B W. *Int J Biochem.* (1989) **21**: 127-137.

Note

This product is for research use only. It is not intended for use in therapeutic or diagnostic procedures for humans or animals. Also, do not use this product as food, cosmetic, or household item, etc. Takara products may not be resold or transferred, modified for resale or transfer, or used to manufacture commercial products without written approval from Takara Bio Inc.

If you require licenses for other use, please contact us by phone at +81 77 565 6973 or from our website at www.takara-bio.com.

Your use of this product is also subject to compliance with any applicable licensing requirements described on the product web page. It is your responsibility to review, understand and adhere to any restrictions imposed by such statements.

All trademarks are the property of their respective owners. Certain trademarks may not be registered in all jurisdictions.

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