For Research Use

TaKaRa

pBApo-EF1α Neo DNA pBApo-EF1α Pur DNA

Product Manual



Table of Contents

| l. | Description | 3 |
|------|------------------------------|---|
| II. | Product Information | 3 |
| III. | Storage | 3 |
| IV. | Vector Map and Cloning Sites | 4 |
| V. | Protocol | 5 |
| VI. | Experimental Examples | 5 |
| VII | Related Products | 6 |

pBApo-EF1α Neo DNA / pBApo-EF1α Pur DNA



I. Description

pBApo-EF1 α is a simple gene expression vector for mammalian cells. This vector carries a promoter from human polypeptide chain elongation factor (EF-1 α promoter) and a polyA signal site from herpes simplex virus thymidine kinase gene. The vectors are useful for the construction of expression plasmids by inserting the ORF of a target gene at the multicloning site. This vector can also be used to express microRNA precursors and other transcripts in addition to ordinary genes.

The pBApo-EF1 α series includes vectors carrying a neomycin-resistance gene or a puromycin-resistance gene as a selection marker in mammalian cells.

II. Product Information

pBApo-EF1 α Neo DNA (Cat. #3243) 20 μ g pBApo-EF1 α Pur DNA (Cat. #3244) 20 μ g

Concentration: $0.5 \mu g/\mu l$

Form: 10 mM Tris-HCl, pH8.0, 1 mM EDTA

III. Storage

-20°C

2 years from date of receipt under proper storage conditions.

IV. Vector Map and Cloning Sites

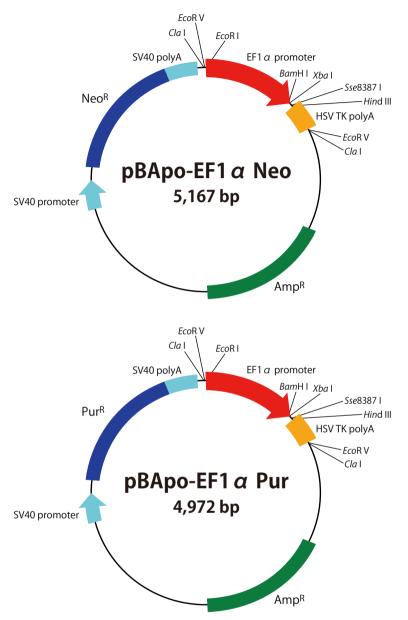


Figure 1. Vector map

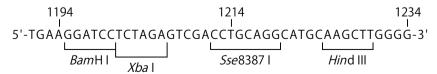


Figure 2. Multicloning site on pBApo-EF1α Neo and pBApo-EF1α Pur

V. Protocol

1. Gene insertion

Insert the ORF of a target gene into the cloning site of a plasmid vector. The ampicillin resistance gene carried by the vector allows for selection of transformed *E. coli*.

2. Transfection

Transfect plasmids using a transfection reagent such as the Xfect[™] series (Cat. #631317 etc.) under the conditions specified in its protocol.

3. Transfected cell selection

- pBApo- EF1 α Neo DNA has a neomycin resistance gene and pBApo- EF1 α Pur DNA has a puromycin resistance gene to allow for drug selection of transfected cells.
- Drug selection should be started at least 24 hours after plasmid transfection. In case
 of a high cell density, reseed cells at appropriate dilution and replace the drug
 containing medium every 3 4 days. Generally, transfected cells can be obtained in
 1 2 weeks.
- Since drug sensitivity varies from cell to cell, determine ahead of time the optimum concentration for the cell used. The concentration will generally be 500 1,000 μ g/ml of G418 for Neo^R gene and 1 3 μ g/ml of puromycin for Pur^R gene.

VI. Experimental Examples

1. Construction of a fluorescent protein expression vector (AcGFP1)

- After digestion of pBApo-EF1α Neo DNA with Hind III, cleaved ends were blunted and further digestion with Xba I was carried out. Subsequently, an approximately 5.1 kb DNA fragment was recovered by agarose gel electrophoresis.
- A DNA fragment of AcGFP1 gene removed from pAcGFP1-C1 Vector (Cat. #632470) by *Nhe* I and *Ssp* I restriction digestions and the digested pBApo-EF1α Neo were ligated using the DNA Ligation Kit <Mighty Mix> (Cat. #6023).
- E. coli JM109 Competent Cells (Cat. #9052) were transformed with the ligation mix and plated on LB plates containing ampicillin.
- The colonies obtained were cultured in 2 5 ml of LB Amp liquid medium to prepare plasmids.
- One of the prepared plasmids was transfected into cultured HeLa cells using Xfect Transfection Reagent (Cat. #631317).
- Cells were observed 24 hours later using a fluorescence microscope to verify the expression of AcGFP1. (Figure 3)

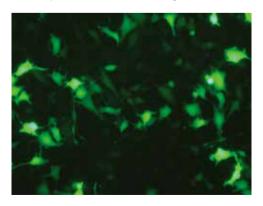


Figure 3. Fluorescence microscopy: 24 hours after transfection of pBApo-EF1a Neo / AcGFP1




2. Comparative expression between EF1a promoter and CMV-IE promoter in mouse ES cells

- Mouse ES E14TG2a cells were transfected with a plasmid (pBApo-EF1α Neo or pBApo-CMV Neo) carrying the AcGFP1 gene using Xfect mESC Transfection Reagent (Cat. #631320).
- Cells were recovered after 48 hours and analyzed for AcGFP1 expression using a flow cytometer (Figure 4. Transient expression).
- Transfected cells were selected by using medium containing 250 μ g/ml G418 and analyzed for AcGFP1 expression using a flow cytometer (Figure 4. Stable expression).

EFp: pBApo-EF1α Neo / AcGFP1 CMVp: pBApo-CMV Neo / AcGFP1 % Positive: the percentage of AcGFP1-positive cells MFI: mean fluorescent intensity of positive cells

Transient expression

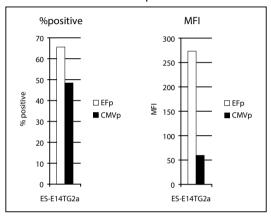

Stable expression

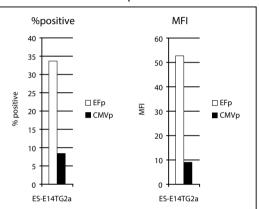


Figure 4. Flow cytometer analyses of transfected cells

Result : The EF1 α promoter provided a higher percentage of AcGFP1-positive cells and a higher level of transient and stable expression in mouse ES cells.

VII. Related Products

pBApo-CMV Vector Series (Cat. #3240_3242)
DNA Ligation Kit < Mighty Mix> (Cat. #6023)

E. coli JM109 Competent Cells (Cat. #9052)

Xfect™ Transfection Reagent (Cat. #631317)

Xfect™ mESC Transfection Reagent (Cat. #631320)
pAcGFP1-C1 Vector (Cat. #632470).
G418 (Cat. #631307, 631308)

Puromycin (Cat. #631305, 631306)



Xfect is a trademark of Takara Bio USA, Inc.

NOTE: This product is for research use only. It is not intended for use in therapeutic or diagnostic procedures for humans or animals. Also, do not use this product as food, cosmetic, or household item, etc.

Takara products may not be resold or transferred, modified for resale or transfer, or used to manufacture commercial products without written approval from TAKARA BIO INC.

If you require licenses for other use, please contact us by phone at +81 77 565 6973 or from our website at www.takara-bio.com.

Your use of this product is also subject to compliance with any applicable licensing requirements described on the product web page. It is your responsibility to review, understand and adhere to any restrictions imposed by such statements.

All trademarks are the property of their respective owners. Certain trademarks may not be registered in all jurisdictions.