pAUR224 DNA

Code No. 3603

Size: 20 µg

Conc.: 1 µg/µl

* 2 years from date of receipt under proper storage conditions.

Description:

pAUR 224 vectors has a mutant *aur1^R* gene derived from *Schizosaccharomyces pombe* as a selective marker that confers Aureobasidin A (AbA)-resistance on cells for yeast transformation system. pAUR224 is a protein expression vector and *E. coli*-yeast shuttle vector. It contains a constitutive promoter of cytomegalovirus (CMV) which function in *Schizo. pombe*.

Form: 10 mM Tris-HCl, pH 8.0
1 mM EDTA

Storage: -20°C

Preparation: cccDNA was purified by CsCl-EtBr ultracentrifugation.

Base pairs: 7,638 bp


Purity:

1. Contains over 60% double-stranded covalently closed circular form I (RFI) DNA.
2. Shown to retain cloning site by dideoxy sequence method.
3. Shown to be cleaved at a single site by restriction enzymes specific to these sites on the cloning site.

Usage:

This is a vector for transformation of yeast *S. pombe* using antibiotic Aureobasidin A. (shuttle vector for protein expression)


Cloning site of pAUR224: bold: stop codon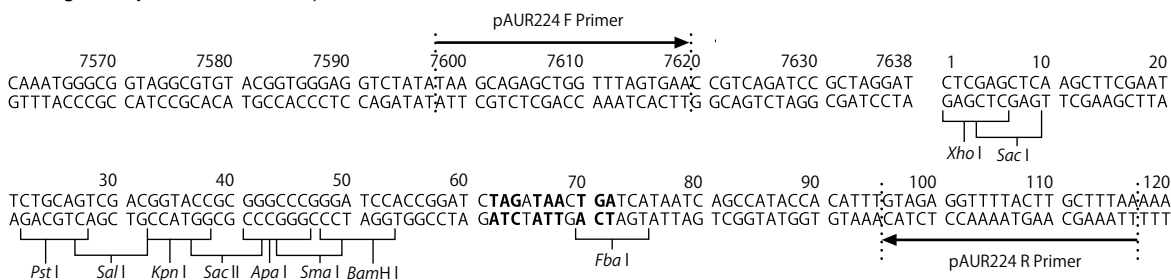

References:

- 1) Takesako K, Kuroda H, Inoue T, Haruna F, Yoshikawa Y, Kato I, Uchida K, Hiratani T, and Yamaguchi H. *J Antibiot.* (1993) **46**: 1414-1420.
- 2) Hashida-Okado T, Yasumoto R, Endo M, Takesako K, and Kato I. *Curr Genet.* (1998) **33**: 38-45.

Vector map of pAUR224 DNA:

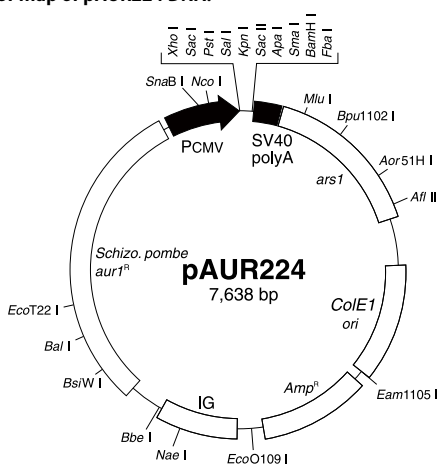- aur1^R*: AbA-resistant gene in *S. pombe*
- ars1*: Replication origin in *S. pombe*
- PCMV: Promoter of CMV
- SV40 polyA: polyA signal from SV40
- Amp^R*: Selective marker in *E. coli*
- ori*: Replication origin in *E. coli*
- IG: intergenic region of M13 phage

Note

This product is for research use only. It is not intended for use in therapeutic or diagnostic procedures for humans or animals. Also, do not use this product as food, cosmetic, or household item, etc. Takara products may not be resold or transferred, modified for resale or transfer, or used to manufacture commercial products without written approval from Takara Bio Inc. If you require licenses for other use, please contact us by phone at +81 77 565 6973 or from our website at www.takara-bio.com. Your use of this product is also subject to compliance with any applicable licensing requirements described on the product web page. It is your responsibility to review, understand and adhere to any restrictions imposed by such statements. All trademarks are the property of their respective owners. Certain trademarks may not be registered in all jurisdictions.

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