# Pfu Protease S

**Code No. 7339**      **Size: 500 U**  
**Conc.: 100 U/ml**

## Description:

*Pfu* Protease S is an endo-type Serine protease with broad specificity toward native and denatured proteins.

## Source:

*Bacillus* species carrying a plasmid that encodes the *Pyrococcus furiosus* protease gene.

**Form:** 25 mM Tris-HCl buffer, pH 7.6 containing 25% ethanol

**Storage:** 4°C

## Quality Control Data :

Please see the Certificate of Analysis (CoA) for each lot. You can download the CoA on Takara Bio website.

**Activity:** 9.5 U/mg (at 95°C)

## Unit definition:

One unit of the enzyme is the amount that produces 1  $\mu$  mol of p-nitroaniline from Suc-AAPF-pNA in a minute at 95°C, pH 7.0.

Suc-AAPF-pNA: N-succinyl Ala-Ala-Pro-Phe p-nitroanilide

## Properties:

Molecular weight: 42,906 Da (amino acid composition)  
45,000 Da (SDS-PAGE)

Optimal temperature: 85 - 95°C

Optimal pH: 6 - 8

Thermal stability: retains 80% activity after 3 hr at 95°C, pH 7.0

## Tolerance to denaturants:

- 1% SDS ; retains 50% activity after 24 hr at 95°C, pH 7.0
- 6.4 M urea ; retains 70% activity after 1 hr at 95°C, pH 7.0
- 50% acetonitrile ; retains 90% activity after 1 hr at 95°C, pH 7.0

Inhibitor: PMSF

Bulk quantities are available. Please contact Takara Bio, Inc. for a bulk quote.

Phone ; +81 77-565-6972      Fax ; +81 77-565-6987

## Note

This product is for research use only. It is not intended for use in therapeutic or diagnostic procedures for humans or animals. Also, do not use this product as food, cosmetic, or household item, etc. Takara products may not be resold or transferred, modified for resale or transfer, or used to manufacture commercial products without written approval from Takara Bio Inc.

If you require licenses for other use, please contact us by phone at +81 77 565 6973 or from our website at [www.takara-bio.com](http://www.takara-bio.com).

Your use of this product is also subject to compliance with any applicable licensing requirements described on the product web page. It is your responsibility to review, understand and adhere to any restrictions imposed by such statements.

All trademarks are the property of their respective owners. Certain trademarks may not be registered in all jurisdictions.

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