# DNA-OFF®

**Code No. 9036**

**Size:**

**500 ml**

## **Description :**

DNA-OFF is a non-alkaline, non-corrosive and noncarcinogenic cleansing solution to eliminate DNA contamination and is intended for use at PCR workstations. Contamination of DNA at a PCR work area can result in DNA amplification artifacts. The solution contains a surfactant. DNA-OFF is stable and heat resistant. DNA-OFF is ready-to-use for eliminating DNA from any surface. By following the simple decontamination instructions below, DNA is completely removed.

## **Storage :**

DNA-OFF should be stored at room temperature. It may generate a precipitate when stored at lower temperatures. The precipitate can be easily brought into solution by incubating at 37°C.

## **Instructions :**

1. Decontamination of laboratory surfaces:  
Apply DNA-OFF directly to the lab surface. Wipe thoroughly with a paper towel, rinse with sterile purified water and dry with a clean paper towel.
2. Decontamination of laboratory apparatus:  
Generously apply DNA-OFF to a paper towel and wipe all exposed surfaces of the apparatus thoroughly. Rinse with a paper towel soaked with sterile purified water and dry with a clean paper towel. To clean small parts, briefly soak them in DNA-OFF, rinse with sterile purified water and dry with a paper towel.
3. Decontamination of plastic and glass vessels:  
Add ample DNA-OFF to coat the entire surface of the vessel by swirling or vortexing. Discard the solution and rinse vessels thoroughly two times with sterile purified water.
4. Decontamination of pipettors:  
Following a manufacturer's instruction, remove the shaft from the pipettor and remove seals and gaskets from the shaft. Soak the shaft for one minute in DNA-OFF, rinse the shaft thoroughly with sterile purified water, let dry and reassemble.  
The other parts than a shaft are cleaned by wiping with a paper towel soaked with DNA-OFF. Then rinse with a paper towel soaked with sterile purified water, and wipe with a clean paper towel.

## **Warning :**

It is recommended to wear plastic gloves when handling DNA-OFF. Contact of solution to skin can cause mild irritation and is harmful if swallowed. In case of contact with eyes, flush with water and contact physician. If swallowed, drink plenty of water and contact physician ; do not induce vomiting.

DNA-OFF is a registered trademark of PureBiotech, LLC.

## **Note**

This product is for research use only. It is not intended for use in therapeutic or diagnostic procedures for humans or animals. Also, do not use this product as food, cosmetic, or household item, etc. Takara products may not be resold or transferred, modified for resale or transfer, or used to manufacture commercial products without written approval from Takara Bio Inc. If you require licenses for other use, please contact us by phone at +81 77 565 6972 or from our website at [www.takarabio.com](http://www.takarabio.com). Your use of this product is also subject to compliance with any applicable licensing requirements described on the product web page. It is your responsibility to review, understand and adhere to any restrictions imposed by such statements. All trademarks are the property of their respective owners. Certain trademarks may not be registered in all jurisdictions.

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