

Takara Bio Europe AB

Cellartis® DEF-CS™ 500 Xeno-Free 3D Spheroid Culture Medium w/o Antibiotics User Manual

Cat. Nos. Y30046, Y30047, Y30048
(061616)

Takara Bio Europe AB
A Takara Bio Company
Arvid Wallgrens backe 20, SE-413 46 Göteborg, Sweden
Europe Technical Support: tech-cellartis@takara-clontech.eu

United States/Canada	Asia Pacific	Europe	Japan
800.662.2566	+1.650.919.7300	+46.(0)31.758.0900	+81.(0)77.543.6116

Page 1 of 9

Table of Contents

I.	Introduction.....	3
II.	List of Components.....	3
III.	Additional Materials Required.....	3
IV.	General Considerations.....	3
	A. Storage and Handling.....	3
	B. Antibiotics.....	4
	C. Cell culture unit.....	4
V.	Culture of Human iPS Cells in Cellartis DEF-CS 500 Xeno-Free 3D Spheroid Culture Medium.....	4
VI.	Cellartis DEF-CS 500 Xeno-Free 3D Spheroid Culture Medium Preparation.....	5
	A. Medium for Passage of Human iPS Cells.....	5
	B. Medium for Maintenance of Human iPS Cells.....	5
VII.	Passage of Human iPS Cells.....	5
	A. Preparations.....	6
	B. Passage.....	6
VIII.	Medium Change for Human iPS Cells.....	6
	A. Preparation.....	6
	B. Medium Change.....	6
IX.	Transfer of Human iPS Cells to Cellartis DEF-CS 500 Xeno-Free 3D Spheroid Culture Medium from other Culture Systems.....	7
	A. Preparation.....	7
	B. Transfer from DEF-CS 500 Xeno-Free Culture Medium.....	7
X.	Cryopreservation of Human iPS Cells.....	8
XI.	Thawing of Human iPS Cells.....	8
XII.	Images of Human iPS Cells Maintained in Cellartis DEF-CS 500 Xeno-Free 3D Spheroid Culture Medium....	8
	Appendix A. Troubleshooting Guide.....	9

Table of Figures

Figure 1.	Schematic presentation of the Cellartis DEF-CS 500 Xeno-Free 3D Spheroid Culture Medium work flow ...	4
Figure 2.	Human iPS cells in Cellartis DEF-CS 500 Xeno-Free 3D Spheroid Culture Medium.....	8
Figure 3.	Human iPS cells in Cellartis DEF-CS 500 Xeno-Free 3D Spheroid Culture Medium (prior to passage, 10X)	8

Table of Tables

Table I.	Guidelines for choice of culture unit (Conical flask size).....	4
Table II.	Guidelines for media volumes at passage and maintenance.....	5
Table III.	Troubleshooting Guide.....	9

I. Introduction

Cellartis DEF-CS 500 Xeno-Free 3D Spheroid Culture Medium is a preclinical grade, chemically defined culture media that is free from human- and animal-derived components and is used for efficient expansion of undifferentiated human induced pluripotent stem (iPS) cells in 3D suspension culture.

All procedures described in the manual have been optimized for Cellartis human iPS cell lines. If you wish to use Cellartis DEF-CS 500 Xeno-Free 3D Spheroid Culture Medium for other human induced pluripotent stem cells, please be aware that procedures and protocols may have to be adjusted.

This product should only be handled by persons who have been trained in laboratory techniques and should only be used in accordance with the principles of good cell culture practice. Takara Bio Europe AB recommends the use of media and reagents according to this manual. Takara Bio Europe AB cannot guarantee correct technical feedback on customer cultures unless the below culture instructions have been followed.

II. List of Components

Cellartis DEF-CS 500 Xeno-Free 3D Spheroid Culture Medium w/o Antibiotics (Cat. No. Y30047)

- **500 ml Cellartis DEF-CS 500 Xeno-Free Basal Medium w/o Antibiotics (Cat. No Y30046; not sold separately)**
- **Cellartis DEF-CS 500 Xeno-Free 3D Spheroid Additives (Cat. No. Y30048)**
 - 500 µl DEF-CS Xeno-Free 3D Spheroid Additive 1 (1,000X)
 - 200 µl DEF-CS Xeno-Free 3D Spheroid Additive 2 (4,000X)
 - 400 µl DEF-CS Xeno-Free 3D Spheroid Additive 3 (500X)

III. Additional Materials Required

The following materials are required but not supplied:

- PBS Dulbecco's w/o Ca²⁺ & Mg²⁺ (D-PBS –/–)
- Accumax (Merck Millipore, Cat. No. SCR006)
- Cell culture vessels, tissue culture treated polystyrene surface, conical flasks or bioreactor
- General cell culture equipment used in cell culture laboratory

If applicable

- Gentamicin (Thermo Fisher, Cat. No. 15750)
- Versene Solution (Thermo Fisher, Cat. No. 15040) (for transfer from monolayer culture)

IV. General Considerations

A. Storage and Handling

Cellartis DEF-CS 500 Xeno-Free Basal Medium w/o Antibiotics (Cat. No. Y30046) should be stored at 4°C; shelf life specified on product label.

Cellartis DEF-CS Xeno-Free 3D Spheroid Additives 1, 2 and 3 should be stored at –20°C; shelf life specified on product label. At first use, thaw provided vials, mix gently and aliquot into appropriate volumes. Store at –20°C according to expiry date on original vial. Thawed vials of Cellartis DEF-CS Xeno-Free 3D Spheroid Additives 1 and 3 may be stored at 4°C for up to one week. Thawed vials of Cellartis DEF-CS Xeno-Free 3D Spheroid Additive 2 should be used the same day. Do not subject the additives to more than a single refreeze and thaw cycle.

NOTE: Cellartis DEF-CS Xeno-Free 3D Spheroid Additive 2 is light sensitive; therefore, avoid unnecessary exposure to light.

B. Antibiotics

If you do elect to use an antibiotic in your medium, we recommend using Gentamicin. Add 0.5 ml to 500 ml of Cellartis DEF-CS 500 Xeno-Free Basal Medium w/o Antibiotics for a final concentration of 50 mg/l.

C. Cell culture unit

This protocol describes how to culture human iPS cell lines in 3D Spheroids in a dynamic system, using a conical flask on an orbital shaker placed in an incubator.

Table I. Guidelines for choice of culture unit (Conical flask size).

Number of seeded cells	Conical flask size	Culture volume
1–4 x 10 ⁶	25 ml	5–15 ml
3–8 x 10 ⁶	100 ml	16–30 ml
7–20 x 10 ⁶	250 ml	35–65 ml

For use of other cell culture vessels or bioreactors, optimization of the protocol will be necessary.

V. Culture of Human iPS Cells in Cellartis DEF-CS 500 Xeno-Free 3D Spheroid Culture Medium

A schematic picture of the transfer, maintenance and cryopreservation of hiPS cell lines in Cellartis DEF-CS 500 Xeno-Free 3D Spheroid Culture Media is shown in Figure 1.

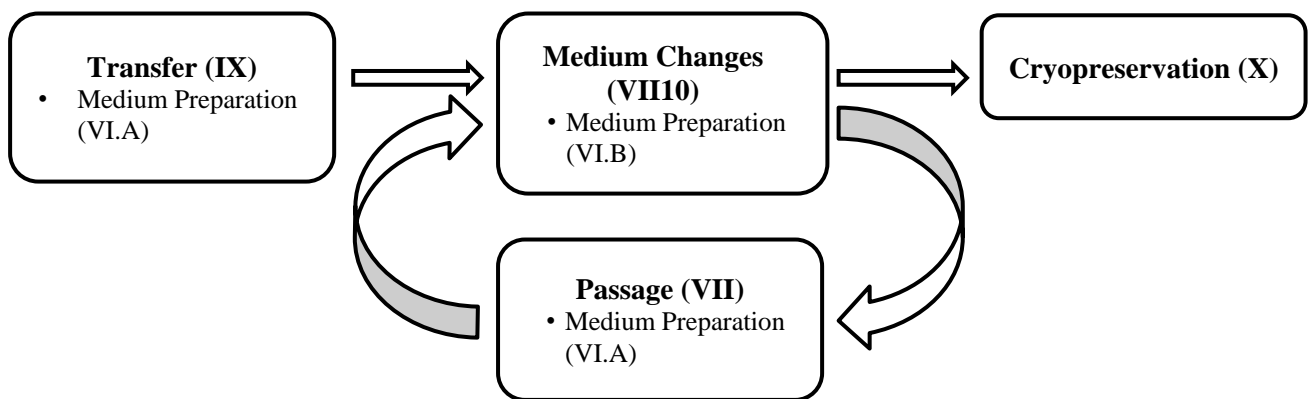


Figure 1. Schematic presentation of the Cellartis DEF-CS 500 Xeno-Free 3D Spheroid Culture Medium work flow. Corresponding sections of this user manual are referenced in brackets.

Human iPS cell lines that are maintained in Cellartis DEF-CS 500 Xeno-Free 3D Spheroid Culture Media should be passaged every 4–5 days for best results. For the first passage from 2D culture into suspension culture, passage at day 7. When the cell density is sparse, you can change the medium every other day; however, it is always important to change medium the day before passage. It is recommended that the cells are grown to a confluence of 1–2 x 10⁶ cells/ml. Do not exceed 3 x 10⁶ cells/ml unless the bioreactor can handle oxygen control and continuous media change (i.e. perfusion system).

VI. Cellartis DEF-CS 500 Xeno-Free 3D Spheroid Culture Medium Preparation

Table II. Guidelines for media volumes at passage and maintenance.

Seeded number of cells	Seeding volume (day 0)	Medium change day 2	Possible medium change day 3	Medium change day before passage	Total volume from seed to harvest	Number of cells at harvest
2×10^6	5–10 ml	10–12 ml	10–12 ml	10–12 ml	36–48 ml	$\sim 10\text{--}15 \times 10^6$
5×10^6	12–25 ml	25–30 ml	25–30 ml	25–30 ml	90–120 ml	$\sim 20\text{--}30 \times 10^6$
10×10^6	25–50 ml	50–60 ml	50–60 ml	50–60 ml	180–240 ml	$\sim 40\text{--}60 \times 10^6$

A. Medium for Passage of Human iPS Cells

1. Decontaminate the external surfaces of all additives and the medium bottle with an appropriate disinfectant and place into the biological safety cabinet.
2. Prepare the appropriate volume of supplemented Cellartis DEF-CS 500 Xeno-Free 3D Spheroid Culture Medium by adding Cellartis DEF-CS Xeno-Free 3D Spheroid Additive 1 (dilute 1:1,000), Additive 2 (dilute 1:4,000) and Additive 3 (dilute 1:500) to prewarmed Cellartis DEF-CS 500 Xeno-Free Basal Medium ($37^\circ\text{C} \pm 1^\circ\text{C}$).
3. Prepare fresh medium on the day of use. Discard any leftover warm medium.

B. Medium for Maintenance of Human iPS Cells

1. Decontaminate the external surfaces of additives and the medium bottle with an appropriate disinfectant and place into the biological safety cabinet.
2. Prepare the appropriate volume of supplemented Cellartis DEF-CS 500 Xeno-Free 3D Spheroid Culture Medium by adding Cellartis DEF-CS Xeno-Free 3D Spheroid Additive 1 (dilute 1:1,000) and Cellartis DEF-CS Xeno-Free 3D Spheroid additive 2 (1:4,000 dilution) to prewarmed Cellartis DEF-CS Xeno-Free Basal Medium ($37^\circ\text{C} \pm 1^\circ\text{C}$). Do not add Cellartis DEF-CS Xeno-Free 3D Spheroid Additive 3 to maintenance medium.
3. Prepare fresh medium on the day of use. Discard any leftover warm medium.

VII. Passage of Human iPS Cells

As a general rule, cells should be seeded at a density of $2\text{--}4 \times 10^5$ cells/ml. Use 2×10^5 cells/ml if leaving the cells five days and $3\text{--}4 \times 10^5$ cells/ml if leaving the cells four days in between passages. However, you should adjust the seeding density to suit your cell line. For recommended culture conditions, see Table I and Table II.

It is recommended that the cells are grown to a confluence of $1\text{--}2 \times 10^6$ cells/ml. If cells are allowed to grow to a higher density, it will have an impact for the next passage, and some cell lines might also be at increased risk of unwanted differentiation. If cultures should appear suboptimal after a few passages, it is recommended to decrease or increase the seeding density. The passage interval may have to be adjusted accordingly.

NOTE: For use of other cell culture vessels or bioreactors, optimization of the protocol will be necessary. As a general rule cells should be seeded at a density of $2\text{--}4 \times 10^5$ cells/ml and grown to a maximum confluence of $2\text{--}3.0 \times 10^6$ cells/ml, unless the bioreactor can handle oxygen control and continuous media change (i.e. perfusion system), in which case higher cell densities can be used.

A. Preparations

Prepare the appropriate volume of supplemented Cellartis DEF-CS 500 Xeno-Free 3D Spheroid Culture Medium for Passage and warm to $37^{\circ}\text{C} \pm 1^{\circ}\text{C}$ as described in Section VI.A. Warm all other reagents to RT.

B. Passage

1. Transfer the spheroid suspension to appropriate centrifuge tubes (15 ml or 50 ml).
2. Centrifuge at 50g for 2 minutes, optimally in a swing-out rotor. Choose a long deceleration time in order to avoid any disturbance in the spheroid pellet due to “forceful braking”.
3. Aspirate the medium and wash the cell spheroids once with D-PBS (–/–). Repeat the centrifugation (step 2). Aspirate the PBS solution.
4. Add 0.1–0.2 ml Accumax/ 10^6 cells (estimate 6-fold increase from seeded number of cells) to the cells and incubate for 15–20 minutes. Tap the side of the tube firmly against the bench 3-5 times to detach cells. If necessary, aspirate the cell suspension 1–3 times with a pipette tip in order to completely disperse the spheroids into a single cell suspension.
5. Dilute the cells in supplemented Cellartis DEF-CS 500 Xeno-Free 3D Spheroid Culture Medium (1:1 dilution).
6. Centrifuge the cells at 200g for 2–3 minutes.
7. Suspend the cells in the appropriate volume of supplemented Cellartis DEF-CS 500 Xeno-Free 3D Spheroid Culture Medium.
8. Count the cells in a haemocytometer or in a cell counter (optimized for the cell type).
9. Add the appropriate volume of cell suspension and medium to the cell culture flask to obtain the selected cell concentration (see Table II).

Note: To minimize the number of spheroids that attach to the walls, pre-wet the conical flask with medium prior to adding the cell suspension.

10. Place the cell culture flask on an orbital shaker at a speed of 60–100 rpm. Place the shaker in an incubator at $37^{\circ}\text{C} \pm 1^{\circ}\text{C}$, 5% CO_2 , and >90% humidity. If necessary, optimize the speed for the specific shaker used.

VIII. Medium Change for Human iPS Cells

When the cell density is sparse, you can change the medium every other day; however, it is always important to change medium the day before passage. If the medium turns yellow due to high metabolic activity, increase the medium volume.

Typically, 3–4 ml medium will be consumed per passage and million cells generated, and as a general assumption cells will expand 4–6 times during each passage. For example, per 1 million cells seeded, you will be able to harvest 4–6 million cells and the total volume of medium consumed for passage and medium changes will be in the range of 18–24 ml. For recommended culture conditions, see Table II.

A. Preparation

Prepare the appropriate volume of supplemented Cellartis DEF-CS 500 Xeno-Free 3D Spheroid Culture Medium for Maintenance as described in Section VI.B and warm to $37^{\circ}\text{C} \pm 1^{\circ}\text{C}$. Warm all other reagents to RT.

B. Medium Change

1. Transfer the spheroid suspension to appropriate centrifuge tubes (15 ml or 50 ml).
2. Separate the spheroids from the bulk of the medium either by using option 1 or option 2.

- *Option 1:* Centrifuge at 50g for 2 minutes, optimally in a swing-out rotor. Choose a long deceleration time in order to avoid any disturbance in the spheroid pellet due to “forceful braking”.
 - *Option 2:* Leave the tube untouched for approximately 5 minutes to allow the spheroids to sink and become concentrated in the bottom of the centrifuge tube.
3. Carefully aspirate the medium.
 4. Suspend the spheroids in prewarmed, supplemented Cellartis DEF-CS 500 Xeno-Free 3D Spheroid Culture Medium and transfer back to the cell culture flask.

Tip: To prevent spheroids from adhering to the plastic surface of the pipette, first aspirate medium up and down one time.

5. Place the cell culture flask on an orbital shaker at a speed of 60–100 rpm. Place the shaker in an incubator at 37°C ± 1°C, 5% CO₂, and >90% humidity. If necessary, optimize the speed for the specific shaker used.

IX. Transfer of Human iPS Cells to Cellartis DEF-CS 500 Xeno-Free 3D Spheroid Culture Medium from other Culture Systems

We recommend transferring your cells to traditional monolayer culture using the Cellartis DEF-CS 500 Xeno-Free Culture Medium (Takara Bio, Cat. No. Y30045), and culturing them in 2D monolayer culture for at least 1–3 passages before transferring to a 3D spheroid culture system.

NOTE: Transferring cells from other culture systems directly into the 3D spheroid culture system is not recommended. However, for some cell lines and culture systems it might be possible to transfer directly into suspension culture. Keep the cell concentration as described for passage (Section VII).

A. Preparation

Prepare the appropriate volume of supplemented Cellartis DEF-CS 500 Xeno-Free 3D Spheroid Culture Medium for Passage, as described in Section VI.A and warm to 37°C ± 1°C. Warm all other reagents to RT.

B. Transfer from DEF-CS 500 Xeno-Free Culture Medium

1. Check cells under microscope; photo document as necessary.
2. Aspirate medium from cell culture vessel and wash the cell layer once with D-PBS (–/–).
3. Add 0.1 ml/cm² of Versene to the cell culture vessel and incubate for 20 minutes or until the cells round up. Detachment should be aided by beating the side of the cell culture vessel firmly or by hitting the short side of the culture vessel against the bench 3–5 times. Avoid flushing the cell layer, if possible.
4. Dilute the cells in supplemented Cellartis DEF-CS 500 Xeno-Free 3D Spheroid Culture Medium (1:1 dilution) and pipette up and down several times to ensure a single cell suspension is achieved.
5. Centrifuge the cells at 300g for 2–5 minutes.
6. Resuspend the cells in the supplemented Cellartis DEF-CS 500 Xeno-Free 3D Spheroid Culture Medium.
7. Count the cells in a haemocytometer or in a cell counter (optimized for the cell type).
8. Add the appropriate volume of cell suspension and medium to the cell culture flask to obtain the selected cell concentration, see Table II.
9. Place the cell culture flask on an orbital shaker at a speed of 60–100 rpm. Place the shaker in an incubator at 37°C ± 1°C, 5% CO₂, and >90% humidity. If necessary, optimize the speed for the specific shaker used.

X. Cryopreservation of Human iPS Cells

Cellartis human iPS cells cultured in Cellartis DEF-CS 500 Xeno-Free 3D Spheroid Culture Medium can be cryopreserved by using common slow freezing protocols for cell suspensions using STEM-CELLBANKER (Takara Bio, Cat. No. CB045). As a general guide, freeze $2.5\text{--}3.5 \times 10^6$ cells in 1 ml freezing medium in a 2 ml cryovial. Freeze the cells as a single cell suspension. Follow steps VII.B 1–9 to achieve a single cell suspension.

NOTE: Cryopreservation of cells in spheroid formation is not recommended.

XI. Thawing of Human iPS Cells

We recommend thawing your cells to traditional monolayer culture using the Cellartis DEF-CS 500 Xeno-Free Culture Medium (Takara Bio, Cat. No. Y30045) and culture them in monolayer culture for at least 1–3 passages before transferring to a 3D Spheroid culture system.

NOTE: Thawing directly into the 3D spheroid culture system is not recommended. However, it may be possible to seed some cell lines directly into suspension culture. Keep the cell concentration as described for passage (see section VII).

XII. Images of Human iPS Cells Maintained in Cellartis DEF-CS 500 Xeno-Free 3D Spheroid Culture Medium

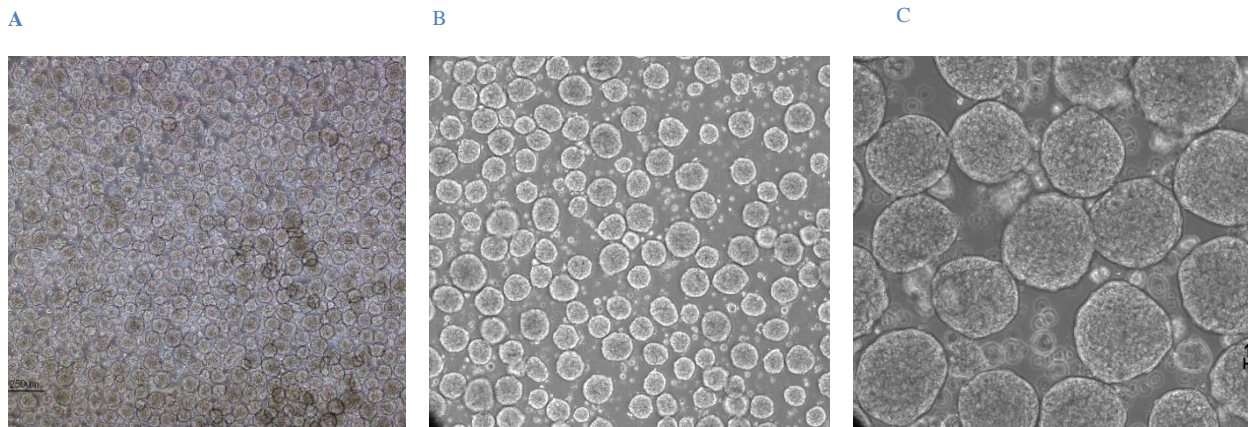


Figure 2. Human iPS cells in Cellartis DEF-CS 500 Xeno-Free 3D Spheroid Culture Medium. Panel A. Day 1 after passage (4X). **Panel B.** Day 1 after passage (10X). **Panel C.** Day 1 after passage (20X).

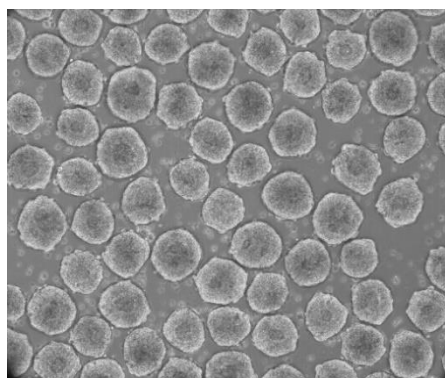


Figure 3. Human iPS cells in Cellartis DEF-CS 500 Xeno-Free 3D Spheroid Culture Medium (prior to passage, 10X).

Appendix A. Troubleshooting Guide

Table III. Troubleshooting Guide

Problem	Possible Explanation	Solution
The cell density at passage varies considerably	Over-compensated cell seeding at previous passages.	Try to keep passage intervals and seeding densities as consistent as possible, i.e. try to not compensate a slow growth for the next passage, or vice versa.
The cells seem to differentiate	Higher concentration of Additive 2 needed.	Increase the additive solution 1.5–2 times compared to the recommended concentration.
The cells seem to differentiate	Too small media volumes used between passages. (Some cell lines have higher metabolic activities, although they do not necessarily divide faster.)	Increase the media volumes used, especially if the medium has turned yellow at higher densities before medium change.
The cells seem to differentiate	Initial cell concentration was too high	Decrease the cell concentration. A good starting concentration is 2×10^5 cells per ml medium, in case the cell line tends to differentiate.
Transferred cells do not adapt to Cellartis DEF-CS 500 Xeno-Free 3D Spheroid Culture Medium	The cells are not used to the new environment.	The cells could benefit from a few passages in Cellartis DEF-CS 500 Xeno-Free Culture Medium prior to transfer to suspension conditions.
Transferred cells do not adapt to Cellartis DEF-CS 500 Xeno-Free 3D Spheroid Culture Medium	The cells are not used to the new environment.	The cells could benefit from a longer initial passage interval. Use a passage interval of 7 days for the first few passages in suspension culture.
The spheres bind to the plastic medium/air interphase	The media volume is not optimal.	It normally helps to increase the media volume 30–50%.
The spheres form big clusters, either directly or after media change	To high initial cell concentration, or non-optimal media volume.	The media volume need to be optimized in conjunction with used shaker platform and cell culture device.

Contact Us	
Customer Service/Ordering	Technical Support
tel: +33.(0)1.3904.6880	tel: +46.(0)31.758.0900
fax: +33.(0)1.3904.6870	fax: +46.(0)31.758.0910
web: www.clontech.com	web: www.clontech.com
e-mail: orders@takara-clontech.eu	e-mail: tech-cellartis@takara-clontech.eu

Notice to Purchaser

This product is for research use only. It is not intended for use in therapeutic or diagnostic procedures for humans or animals. Also, do not use this product as food, cosmetic, or household item, etc.

This product may not be resold or transferred, modified for resale or transfer, or used to manufacture commercial products without written approval from Takara Bio Europe AB.

If you require licenses for other use, please contact us by phone at +46 31 758 0900.

Your use of this product is also subject to compliance with any applicable licensing requirements as detailed in our catalogues, on our website at <http://www.clontech.com>, on the label or other documentation accompanying the goods. It is your responsibility to review, understand and adhere to any restrictions imposed by such statements.

Takara and the Takara logo are trademarks of TAKARA HOLDINGS INC., Kyoto, Japan. Cellartis and DEF-CS are trademarks of Takara Bio Europe AB. All other trademarks are the property of their respective owners. Certain trademarks may not be registered in all jurisdictions. Takara Bio Europe AB is a Takara Bio company. ©2016 Takara Bio Europe AB

This document has been reviewed and approved by the Takara Bio Europe AB Quality Assurance Department.