Yatalase

Code No. T017

Size : 2 g

Example of protoplast preparation :

- (1) *Aspergillus oryzae*

| | |
|------------|---|
| Medium | Czapek-Dox + 0.5% Casamino acid (pH 5.6) |
| Culture | 30°C, 20 hr, Shake culture (Rotary at 140 rpm) |
| Conditions | 2% Yatalase solution 0.6 M (NH ₄) ₂ SO ₄ 50 mM Maleate Buffer (pH 5.5) 30°C Reciprocal shaking (at 60 - 70 rpm) 2 - 3 hr |
- (2) *Aspergillus kawachi*

| | |
|------------|---------------------------------|
| Medium | Same as in (1) |
| Culture | 30°C, 20 hr, Stationary culture |
| Conditions | Same as in (1) |
- (3) *Aspergillus terreus, Penicillium citrinum, Penicillium lanosum, Tricoderma koningii*

| | |
|------------|--------------------------|
| Medium | Dextrin-peptone (pH 5.5) |
| Culture | Same as in (1) |
| Conditions | Same as in (1) |
- (4) *Mucor hiemalis, Rhizopus nigricans*

| | |
|------------|---|
| Medium | 2% Malt extract |
| Culture | 30°C, 12 hr, Stationary culture (Sporangiospores in germ) |
| Conditions | 2% Yatalase solution 0.5 M MgSO ₄ 50 mM Maleate Buffer (pH 5.5) 30°C Reciprocal shaking (at 60 - 70 rpm) 4 hr |
- (5) *Pleurotus ostreatus, Coprinus cinereus, Lentinus edodes*

| | |
|------------|---|
| Medium | OSG or MYG medium (pH 5.5) |
| Culture | 25 - 30°C, 3 - 4 days, Stationary culture |
| Conditions | 2% Yatalase solution 0.6 M MgSO ₄ 50 mM Maleate Buffer (pH 5.5) 30°C Reciprocal shaking (at 60 - 70 rpm) 2 - 3 hr |
- (6) *Monascus sp.*

| | |
|------------|--|
| Medium | Same as in (3) |
| Culture | 25°C, 20 hr, Shake culture (Rotary at 140 rpm) |
| Conditions | Same as in (1) |

Description :

This product was prepared from lipoid culture supernatant of *Corynebacterium* sp. OZ-21. This product has complex lytic activities of fungal cell mainly consisting of chitinase and chitobiase activity.

Origin : *Corynebacterium* sp. OZ-21

Appearance : Lyophilized powder (containing lactose as the excipient)

Storage : 4°C, dry condition

Unit definition :

[Chitinase activity]

1 unit is defined as the amount of enzyme required to release 1 μ mol of *N*-Acetylglucosamine from 0.5% chitin powder solution in 1 min at 37°C, pH 6.0.

[Chitobiase activity]

1 unit is defined as the amount of enzyme required to release 1 μ mol of *p*-Nitrophenol from *p*-Nitrophenyl-*N*-acetyl- β -D-glucosaminide solution in 1 min at 37°C, pH 6.0.

[Lytic activity]

1 unit is defined as the amount of enzyme required to cause a 1% decrease in absorbance at 660 nm in 1 hour under the assay conditions described below.

Solutions

Solution A (Substrate solution) ;

Harvest a culture of *Aspergillus oryzae* (cultured in dextrin-peptone medium at 25 to 30°C for 1 to 2 days) and homogenize on a Waring blender (16,000 rpm, 3 min). Subsequently, apply a French press process to produce fragments. Then, wash the fragments thoroughly with water over filter paper, and then with acetone, followed by diethyl ether. Air-dry the pellet and suspend in 0.1 M acetate buffer, pH 6.0, to give 0.125% (w/v).

Solution B (Enzyme solution);

Dissolve the enzyme in sterile purified water (0.8 mg/ml).

Assay system

1. Transfer 4 ml of solution A to a test tube.
2. Add 1 ml of solution B and immediately vortex.
3. Incubate at 37°C by gently shaking for 1 hr.
4. Read the absorbance at 660 nm.

Quality Control Data :

Please see the Certificate of Analysis (CoA) for each lot. You can download the CoA on Takara Bio website.

Manufactured by Ozeki corporation.

Note

This product is for research use only. It is not intended for use in therapeutic or diagnostic procedures for humans or animals. Also, do not use this product as food, cosmetic, or household item, etc. Takara products may not be resold or transferred, modified for resale or transfer, or used to manufacture commercial products without written approval from Takara Bio Inc. If you require licenses for other use, please contact us by phone at +81 77 565 6973 or from our website at www.takara-bio.com. Your use of this product is also subject to compliance with any applicable licensing requirements described on the product web page. It is your responsibility to review, understand and adhere to any restrictions imposed by such statements. All trademarks are the property of their respective owners. Certain trademarks may not be registered in all jurisdictions.

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