

pOET1C 6xHis Transfer Plasmid

Product Data Sheet

Product information for MIR 6151

SDS and Certificate of Analysis available at mirusbio.com/literature



SPECIFICATIONS

Storage	Store pOET1C_6xHis Transfer Plasmid DNA at -20°C.
Product Guarantee	1 year from the date of purchase, when properly stored and handled.
Concentration	20 µl plasmid DNA at 500 ng/µl, sterile filtered in TE Buffer.

Plasmid Description: pOET1C_6xHis, officially pOET1.1C_6xHis, is a baculovirus transfer vector designed to facilitate high level expression of foreign genes under the AcMNPV polyhedrin promoter in insect cells. Following transfection into insect cells, the AcMNPV sequences flanking the expression cassette allow recombination into the baculovirus genome which replicates to produce baculovirus that can be harvested directly from the cell culture medium. pOET1C_6xHis plasmid features include a Col E1 origin of replication, an ampicillin resistance gene for selection in *E. coli*, a multiple cloning site (MCS) for insertion of the foreign gene in the correct frame and orientation and a C-terminal 6xHis-Tag[®] fusion sequence that may be utilized for recombinant protein purification with Ni-NTA agarose columns and removed by subsequent incubation with the proteolytic enzyme, Thrombin. pOET1C_6xHis is compatible with the *flashBAC*[™] System or any baculovirus expression system that utilizes homologous recombination at the polyhedrin locus in insect cells.

Plasmid Features:

AcMNPV: 623 - 1769

Polyhedrin (PH) promoter: 1770 - 1870


Multiple Cloning Site (MCS): 1908 - 1982

6xHis: 1982 - 2000

AcMNPV: 2012 - 2628

Col E1 Replication Origin: 2834 - 3453

Amp Resistance Marker: 3608 - 4468

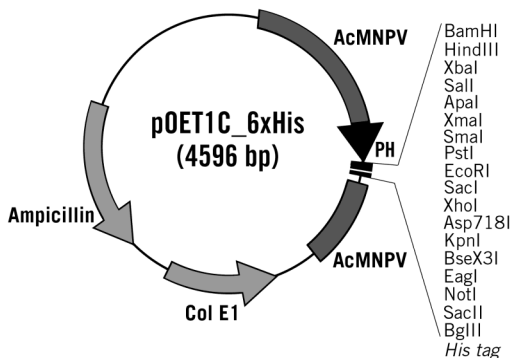

Additional Information: All pOET vectors can be propagated in DH5α or other general purpose *E. coli* strains. For more information on baculovirus expression using the *flashBAC*[™] System, pOET transfer plasmids or optimized insect cell transfections using *TransIT*[®]-Insect Transfection Reagent, please visit www.mirusbio.com.

For Research Use Only

Mirus Bio LLC

www.mirusbio.com | techsupport@mirusbio.com | U.S. Toll Free: 844.647.8724 | Direct: +1.608.441.2852

► NOTES



SDS and Certificate of Analysis available at mirusbio.com/literature

flashBAC[™] Systems and pOET vectors are sold by Mirus Bio through partnership with Oxford Expression Technologies, Oxford, UK.

©1996-2024 All rights reserved. Mirus Bio LLC. All trademarks are the property of their respective owners.
For terms and conditions, visit www.mirusbio.com

Rev2 13SEP2024

Mirus Bio LLC

www.mirusbio.com | techsupport@mirusbio.com | U.S. Toll Free: 844.647.8724 | Direct: +1.608.441.2852