

RevIT™ AAV Enhancer

Quick Reference Protocol

Instructions for MIR 8000, 8006

SDS and Certificate of Analysis available at mirusbio.com/8000



SPECIFICATIONS

Storage	Store RevIT™ AAV Enhancer at -10 to -30°C, tightly capped. Before each use , warm to room temperature and vortex gently. Do not freeze/thaw > 5 times.
Product Guarantee	When properly stored and handled, RevIT™ AAV Enhancer is guaranteed for 6 months from the date of purchase.

RevIT™ AAV Enhancer Workflow

Maintain Cells

Passage cells regularly and ensure they are >95% viable before transfection.

Day 0

Seed cells.

Per ml of culture, add **RevIT™**: 0.5 - 1.5 µl to diluted DNA prior to adding the transfection reagent as directed by the manufacturer.

Day 2 - 3

Harvest AAV **48 - 72 hr** post-transfection.

Product Description

Recombinant AAV has become an invaluable tool for gene therapy and the creation of isogenic human disease models. RevIT™ AAV Enhancer boosts the performance of transfection reagents used in AAV production applications or workflows. The RevIT™ AAV Enhancer is ideal for generating high titer AAV preparations to accelerate research and development.

Product Formats

RevIT™ AAV Enhancer is supplied in the following formats:

Product No.	Quantity
MIR 8000	1.5 ml
MIR 8006	10 × 1.5 ml

RevIT™ AAV Enhancer can be paired with TransIT-VirusGEN® Transfection Reagent in 3 ml (MIR 8007) and 30 ml (MIR 8008) configurations.

Product Usage

Mix RevIT™ AAV Enhancer well prior to use. Perform transfection per transfection reagent manufacturer's recommendations. Prior to adding transfection reagent, add RevIT™ AAV Enhancer to diluted DNA.

Titrate RevIT™ AAV Enhancer from 0.5 to 1.5 µl per 1 ml of cell culture media to determine the optimal amount for production of your specific viral vector. For example, if transfecting a 30 ml cell culture, test 15 µl to 45 µl of RevIT™ AAV Enhancer.

Harvest AAV as appropriate for your process, typically 48 - 72 hours after transfection.

In need of higher titers and percent full capsids?

Pair RevIT™ AAV Enhancer with TransIT-VirusGEN® to yield even better performance.

Not sure? Request a free sample of TransIT-VirusGEN® here: mirusbio.com/sample

For Research Use Only

Mirus Bio LLC

www.mirusbio.com | techsupport@mirusbio.com | U.S. Toll Free: 844.647.8724 | Direct: +1.608.441.2852



Reagent Agent[®]

Reagent Agent[®] is an online tool designed to help determine the best solution for nucleic acid delivery based on in-house data, customer feedback and citations.

Learn more at: [mirusbio.com/ra](https://www.mirusbio.com/ra)

©1996-2023 All rights reserved. Mirus Bio LLC. All trademarks are the property of their respective owners.
For terms and conditions, visit www.mirusbio.com

Rev0 070723

Mirus Bio LLC

www.mirusbio.com | techsupport@mirusbio.com | U.S. Toll Free: 844.647.8724 | Direct: +1.608.441.2852