50 bp DNA Ladder (Dye Plus)

Code No. 3421A **Size:** 500 µl
 (for 100 lanes)
Conc.: 200 ng/µl

Description:

This marker consists of 10 fragments between 50 and 500 bp in multiples of 50 bp and additional fragments at 600, 700, 800, 900, 1,000, and 1,500 bp. The 300 and 600 bp fragments are brighter than the other fragments and serve as a visible reference indicator; all other fragments appear with equal intensity on the gel. They are all double-stranded DNA fragments. Since this product is premixed with loading dyes and glycerol, it is possible to apply this product to agarose gel electrophoresis directly.

Form:

8.33 mM	Tris-HCl, pH 8.0
5.83 mM	EDTA
6%	Glycerol
0.0083%	Bromophenol Blue
0.0033%	Xylene Cyanol

Storage:

-20°C
Once thawed, this product should be stored at 4°C.
Avoid freeze-thaw cycles.

Note: Once opened or thawed, store at 4°C and use earlier avoiding contamination.

Usage:

Usage:
Used as a DNA molecular size marker in agarose gel electrophoresis.

Application example:

Apply 5 μ l of DNA Ladder on agarose gel. Run agarose gel electrophoresis using ex. 3% PrimeGel™ Agarose PCR-Sieve (Cat. #5810A). Perform staining with EtBr, or SYBR™ Green I Nucleic Acid Gel Stain (Cat. #5760A/5761A)*.

* Not available in all geographic locations. Check for availability in your area.

Notice:

Notice:
This product is designed for the use with agarose gel.
When the marker is used for other purpose, the separation pattern may be strange. When marker is used for denatured polyacrylamide gel electrophoresis, duplicated bands may appear in a gel and accurate isolation cannot be performed. When used for non-denatured polyacrylamide gel, this product may also show duplicated bands under some condition. Avoid using this product for polyacrylamide gel electrophoresis.

PrimeGel is a trademark of Takara Bio Inc.
SYBR is a trademark of Life Technologies Corporation.

Note

NOTE This product is for research use only. It is not intended for use in therapeutic or diagnostic procedures for humans or animals. Also, do not use this product as food, cosmetic, or household item, etc. Takara products may not be resold or transferred, modified for resale or transfer, or used to manufacture commercial products without written approval from Takara Bio Inc. If you require licenses for other use, please contact us from our website at www.takarabio.com. Your use of this product is also subject to compliance with any applicable licensing requirements described on the product web page. It is your responsibility to review, understand and adhere to any restrictions imposed by such statements. All trademarks are the property of their respective owners. Certain trademarks may not be registered in all jurisdictions.

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