

Cat. # 9028

For Research Use

TAKARA

***E. coli* HST08 Premium
Electro-Cells**

Product Manual

v202012Da

Table of Contents

I.	Description.....	3
II.	Components	3
III.	Storage	3
IV.	Protocol.....	4
V.	Quality.....	6
VI.	Genotype.....	6
VII.	Cell Density	6
VIII.	References.....	6
IX.	Related Products	6

I. Description

E. coli HST08 Premium Electro-Cells are specially prepared to be highly appropriate to electroporation. Electroporation is used to transfer DNA into cells by perforating the cytoplasmic membrane with a high voltage pulse.

In addition, *E. coli* HST08 Premium Electro-Cells lack the genes necessary for digesting foreign methylated DNA: *mrr*, *hsdRMS*, *mcrBC*, and *mcrA*. These properties make these competent cells useful in a wide range of applications such as methylated DNA cloning, genetic library construction, and BAC-based long-length genome libraries. Even when working with very large plasmids, the transformation efficiency and colony growth rates remain very high*. Effective DNA cloning and genetic library construction with insert DNA larger than 10 kb is possible if this product is used with TaKaRa DNA Ligation Kit LONG (Cat. #6024).

* Compared with other competent cells of the same genotype.

For transformation of pUC-based plasmids, selection of recombinants may be simplified by adding X-Gal to the media, thus using the α -complementarity to β -galactosidase of the competent cell to visualize vector-insert containing transformants.

X-Gal : 5-Bromo-4-Chloro-3-Indolyl- β -D-Galactoside

II. Components

<i>E. coli</i> HST08 Premium Electro-Cells	50 μ l x 10
pUC19 plasmid (0.1 ng/ μ l)	10 μ l
SOC Medium*	1 ml x 10

* SOC Medium :	2%	Tryptone
	0.5%	Yeast extract
	10 mM	NaCl
	2.5 mM	KCl
	10 mM	MgSO ₄
	10 mM	MgCl ₂
	20 mM	Glucose

III. Storage

-80°C

Note : Store at -80°C or lower. If the storage temperature is not maintained consistently, the transformation efficiency will be reduced. You may determine the transformation efficiency of stored cells by using the included pUC19 control.

Do not store in liquid nitrogen.

IV. Protocol

- (1) Thaw 50 μ l of *E. coli* HST08 Premium Electro-Cells on ice just before use.
 - (2) Add 1 - 2 μ l of DNA solution*¹ into the thawed cell suspension.
 - (3) Transfer the mixture of cells and DNA to a cold 0.1 cm electroporation cuvette.
 - (4) After applying a pulse*², immediately add 1 ml of SOC Medium (prechilled on ice).
 - (5) Incubate at 37°C for 1 hour with shaking at 160 - 225 rpm.
 - (6) Plate an appropriate amount of culture*³.
 - (7) Place plates in a 37°C incubator and grow overnight.
- * 1 When the sample DNA solution contains salt(s), dilute it with TE buffer or sterile purified water. Alternatively, desalination by ethanol precipitation is recommended. Ethanol precipitation should be performed when the sample DNA is prepared for use with the TaKaRa DNA Ligation Kit.
 - * 2 Takara Bio uses BIO-RAD MicroPulser and the electrical condition is 1.5 kV. In the case of BIO-RAD Gene Pulser, standard electrical conditions are 200 Ω , 25 μ F and 1.5 kV.
 - * 3 Plate no more than 100 μ l for a ϕ 9 cm plate. If necessary, dilute the culture with the same medium as used in step (4).

[Please read before proceeding]

1. Place a tube of Electro-Cells in a dry ice / EtOH bath immediately upon removal from the -80°C freezer. Keep the cells in the bath until you are ready to proceed.
2. For 50 μ l of Electro-Cells, use no more than 10 ng of high purity DNA, or the transformation efficiency may be reduced.
3. For 50 μ l of Electro-Cells, use no more than 5 μ l of DNA solution, or the transformation efficiency may be reduced.
4. If you change the quantity of competent cells electroporated, it may be necessary to reevaluate the conditions.
5. L-broth or ϕ b-broth can be used instead of SOC Medium, but efficiency may be reduced.

<u>L-broth</u> :	<u>Ingredient</u>	<u>per liter water</u>
	Tryptone	10 g
	Yeast extract	5 g
	NaCl	5 g

Adjust to around pH 7.5 with 1 N NaOH and autoclave.

<u>ϕb-broth</u> :	<u>Ingredient</u>	<u>per liter water</u>
	Tryptone	20 g
	Yeast extract	5 g
	MgSO ₄ ·7H ₂ O	5 g

Adjust to around pH 7.5 with 1 N KOH and autoclave.

6. When using X-Gal, follow the procedures described as below:
Add 20 mg/ml X-Gal (dissolved in dimethylformamide) at a ratio of 200 - 300 μ l/100 ml agar media.
7. Once the competent cells have been thawed, refreezing for storage is not recommended. If this is unavoidable, flash freeze the cells on dry ice/ethanol and store them promptly at -80°C. However, the transformation efficiency will be lowered by at least one order of magnitude.

V. Quality

1. Efficiency of transformation
10 pg of pUC19 plasmid was transformed according to section IV and transformants were selected on a L-plate containing ampicillin. The resulting transformation efficiency was $> 1 \times 10^9$ colonies / μ g pUC19 plasmid.
2. Confirmation of β -galactosidase α -complementation.
When transformed with a pUC19 plasmid, blue colonies appeared on a L-agar plate containing 100 μ g/ml ampicillin, and 60 μ g/ml X-Gal.

VI. Genotype

E. coli HST08 Premium :
F⁻, *endA1*, *supE44*, *thi-1*, *recA1*, *relA1*, *gyrA96*, *phoA*, Φ 80d *lacZ* Δ M15,
 Δ (*lacZYA -argF*) U169, Δ (*mrr - hsdRMS - mcrBC*), Δ *mcrA*, λ ⁻

VII. Cell density

$> 1 \times 10^{10}$ bacteria/ml

VIII. References

- 1) Dower W J, Miller J F, and Ragsdale C W. *Nucl Acids Res.* (1988) **16**: 6127.
- 2) Bottger E C. *Biotechniques.* (1988) **6**: 878.

IX. Related Products

E. coli HST08 Premium Competent Cells (Cat. #9128)
TaKaRa DNA Ligation Kit LONG (Cat. #6024)
X-Gal (5-Bromo-4-Chloro-3-Indolyl- β -D-Galactoside) (Cat. #9031)
pUC19 DNA (Cat. #3219)

NOTE: This product is for research use only. It is not intended for use in therapeutic or diagnostic procedures for humans or animals. Also, do not use this product as food, cosmetic, or household item, etc.

Takara products may not be resold or transferred, modified for resale or transfer, or used to manufacture commercial products without written approval from Takara Bio Inc.

If you require licenses for other use, please contact us by phone at +81 77 565 6972 or from our website at www.takara-bio.com.

Your use of this product is also subject to compliance with any applicable licensing requirements described on the product web page. It is your responsibility to review, understand and adhere to any restrictions imposed by such statements.

All trademarks are the property of their respective owners. Certain trademarks may not be registered in all jurisdictions.
