Anti-Human Osteocalcin, Monoclonal (Clone 5-12H)

Code No. M184

Size : 0.1 mg Mouse Ig

Subclass : IgG2b

* 2 years from date of receipt under proper storage conditions.

Source :

Monoclonal antibody was obtained by fusing the mouse myeloma cell-line P3U1 with spleen cells of C57BL/6 mouse after immunization with KLH-conjugated peptide from human osteocalcin (amino acids 1-6 YLYQWL). The monoclonal antibody was harvested from ascitic fluid of SCID mouse.

Purification :

Antibody was purified by column chromatography, dissolved in 10 mM PBS, pH 7.4, containing 1.0% bovine serum albumin, and then lyophilized. The lyophilized antibody does not contain preservative.

Form : Lyophilized

Reconstitution :

Dissolve the lyophilized antibody in 50 μ l of distilled water (final concentration: 2.0 mg/ml). This solution can be used as a stock solution. If dilution is necessary for your application, dilute the stock solution with the following Dilution solution just prior to use. When the entire amount of antibody is to be used over a short time period, it may be dissolved directly in 500 μ l or more of the Dilution solution.

Note (1) : Be sure to store the antibody at a minimum concentration of 2.0 mg/ml. A lower antibody concentration may result in decreased stability.

Note (2) : Reconstituted antibody solution should contain 0.1% sodium azide as a preservative when stored at 4°C.

Dilution solution:

10 mM PBS (pH 7.4)
1.0% BSA
(0.1% NaN_3) *

* When stored at 4°C, 0.1% sodium azide should be added as a preservative.

Specificity:

- This antibody recognizes amino acid sequence (1 - 6) of human osteocalcin.
- This antibody reacts with the full sequence of human osteocalcin.

Cross reactivity :

This antibody does not react with bovine, rat, guinea pig, pig, horse, pigeon and goose osteocalcin.

Storage : 4°C

This product does not contain preservative.

The stock solution (2.0 mg/ml) should be stored in aliquots at -20°C for 1 year, or should be stored at 4°C for 6 months after adding 0.1% sodium azide. Avoid repeated freeze-thaw cycles. Diluted antibody should not be stored.

Working concentration:

- 1 - 10 μ g/ml (for Western blotting by color detection)
- 20 - 40 μ g/ml (for immunohistochemical staining by color detection)

Application:

- Immunohistochemical detection of osteocalcin on formalin fixed, decalcified, paraffin embedded human tissue sections
 - * Antigen retrieval: Proteinase K treatment
- Western blot analysis of human osteocalcin under reducing or non-reducing conditions

Additional information :

Sandwich immunoassay is possible by combination of clones 5-12H (solid phase) and GlaOC4-30 (labelled). However this product is not suitable for Sandwich ELISA assay, because it contains BSA.

Note

This product is for research use only. It is not intended for use in therapeutic or diagnostic procedures for humans or animals. Also, do not use this product as food, cosmetic, or household item, etc. Takara products may not be resold or transferred, modified for resale or transfer, or used to manufacture commercial products without written approval from Takara Bio Inc. If you require licenses for other use, please contact us by phone at +81 77 565 6973 or from our website at www.takara-bio.com. Your use of this product is also subject to compliance with any applicable licensing requirements described on the product web page. It is your responsibility to review, understand and adhere to any restrictions imposed by such statements. All trademarks are the property of their respective owners. Certain trademarks may not be registered in all jurisdictions.

Anti-Human Osteocalcin, Monoclonal (Clone 5-12H)

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