Anti-ProS2, Monoclonal

Code No. M200 **Size : 0.1 mg Mouse Ig**
Clone No. ProS 7B-8F **Subclass : IgG1**

* 2 years from date of receipt under proper storage conditions.

Source:

Monoclonal antibody was obtained by fusing the mouse myeloma cell-line P3U1 with lympho cell of C57BL/6 mouse after immunization with ProS2 tag protein. The monoclonal antibody was harvested from ascitic fluid of SCID mouse.

* ProS2 is solubilization tag of about 23 kDa, which is tandem-dimer of N-terminal domain of a soluble protein, Protein S, derived from myxobacteria *Myxococcus xanthus*. It is used at *E. coli* cold shock expression system pCold™ ProS2 DNA (Cat. #3371).

Purification:

Antibody was purified by column chromatography, dissolved in 10 mM PBS, pH 7.4, containing 1.0% bovine serum albumin, and then lyophilized. The lyophilized antibody does not contain preservative.

Form: Lyophilized

Reconstitution:

Dissolve the lyophilized antibody in 50 μ l of distilled water (final concentration 2.0 mg/ml). This solution can be used as a stock solution. If dilution is necessary for your application, dilute the stock solution with the following Dilution solution just prior to use. When the entire amount of antibody is to be used over a short time period, it may be dissolved directly in 500 μ l or more of the Dilution solution.

Note (1) : Be sure to store the antibody at a minimum concentration of 2.0 mg/ml. A lower antibody concentration may result in decreased stability.

Note (2) : Reconstituted antibody solution should contain 0.1% sodium azide as a preservative when stored at 4°C.

Dilution solution:

10 mM PBS (pH 7.4)
1.0% BSA
(0.1% NaN₃)*

* When stored at 4°C, 0.1% sodium azide should be added as a preservative.

Specificity:

This antibody specifically reacts with ProS2 tag of a fusion protein expressed with pCold ProS2 DNA.

Cross reactivity:

This antibody does not react with proteins derived from *E. coli*, and TF (Trigger Factor) tag.

Storage: 4°C

This product does not contain preservative.

The stock solution (2.0 mg/ml) should be stored in aliquots at -20°C for 1 year, or should be stored at 4°C for 6 months after adding 0.1% sodium azide. Avoid repeated freeze-thaw cycles. Diluted antibody should not be stored.

Working concentration :

- Western Blotting by color detection : 0.5 - 1 μ g/ml
- Immunoprecipitation : 10 μ g/ml

Application:

For detection of ProS2-fused protein expressed by using pCold ProS2 DNA

- Western blot analysis under reducing or non-reducing conditions
- Immunoprecipitation

References:

Significant Enhanced Expression and Solubility of Human Proteins in *E. coli* by Fusion with Protein S from *Myxococcus xanthus*. Kobayashi H, Yoshida T, and Inouye M *Appl Environ Microbiol.* (2009) **75**: 5356-5362.

pCold is a trademark of Takara Bio Inc.

Note

This product is for research use only. It is not intended for use in therapeutic or diagnostic procedures for humans or animals. Also, do not use this product as food, cosmetic, or household item, etc. Takara products may not be resold or transferred, modified for resale or transfer, or used to manufacture commercial products without written approval from Takara Bio Inc. If you require licenses for other use, please contact us by phone at +81 77 565 6973 or from our website at www.takara-bio.com. Your use of this product is also subject to compliance with any applicable licensing requirements described on the product web page. It is your responsibility to review, understand and adhere to any restrictions imposed by such statements. All trademarks are the property of their respective owners. Certain trademarks may not be registered in all jurisdictions.

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