Anti-TF, Monoclonal

Code No. M201 **Size : 0.1 mg Mouse Ig**
Clone No. TF 19-7F **Subclass : IgG1**

* 2 years from date of receipt under proper storage conditions.

Source:

Monoclonal antibody was obtained by fusing the P3U1 myeloma cell-line with lymph node cells of C57BL/6 mouse after immunization with synthetic peptide (165-178) [TIDFTGSVDGEEFE] of Trigger Factor * conjugated with KLH. The monoclonal antibody was harvested from ascitic fluid of SCID mouse.

* Trigger Factor (TF) is a chaperon protein derived from *E. coli*. It is used at *E. coli* cold shock expression system pCold™ TF DNA (Cat. #3365).

Gene : tig

Protein : peptidyl-prolyl cis/trans isomerase (trigger factor)

Origin : [*Escherichia coli* str. K-12 substr. MG1655]

Purification :

Antibody was purified by column chromatography, dissolved in 10 mM PBS, pH 7.4, containing 1.0% bovine serum albumin, and then lyophilized.

The lyophilized antibody does not contain preservative.

Form: Lyophilized

Reconstitution:

Dissolve the lyophilized antibody in 50 μ l of distilled water (final concentration 2.0 mg/ml). This solution can be used as a stock solution. If dilution is necessary for your application, dilute the stock solution with the following Dilution solution just prior to use. When the entire amount of antibody is to be used over a short time period, it may be dissolved directly in 500 μ l or more of the Dilution solution.

Note (1) : Be sure to store the antibody at a minimum concentration of 2.0 mg/ml. A lower antibody concentration may result in decreased stability.

Note (2) : Reconstituted antibody solution should contain 0.1% sodium azide as a preservative when stored at 4°C.

Dilution solution:

10 mM PBS (pH 7.4)

1.0% BSA

(0.1% NaN₃)*

* When stored at 4°C, 0.1% sodium azide should be added as a preservative.

Specificity:

This antibody specifically reacts with Trigger Factor (TF) tag of a fusion protein expressed with pCold TF DNA.

Cross reactivity :

This antibody does not react with other protein derived from *E. coli*, and ProS2 tag.

Storage: 4°C

This product does not contain preservative.

The stock solution (2.0 mg/ml) should be stored in aliquots at -20°C for 1 year, or should be stored at 4°C for 6 months after adding 0.1% sodium azide. Avoid repeated freeze-thaw cycles. Diluted antibody should not be stored.

Working concentration :

- Western Blotting by color detection : 1 - 5 μ g/ml
- Immunoprecipitation : 5 - 10 μ g/ml *

* Attach 5 - 10 μ g of this antibody to 50 - 100 μ l of anti-Mouse IgG resin (or beads). Mix the 1 ml of sample and the immobilized antibody resin (beads) well. Wash the immune-complex resin (beads) and analyze by SDS-PAGE or Western Blotting.

Application :

- For detection of TF-fused protein expressed by using pCold TF DNA
- Western blot analysis under reducing and non-reducing conditions.
 - Immunoprecipitation

pCold is a trademark of Takara Bio Inc.

Note

This product is for research use only. It is not intended for use in therapeutic or diagnostic procedures for humans or animals. Also, do not use this product as food, cosmetic, or household item, etc. Takara products may not be resold or transferred, modified for resale or transfer, or used to manufacture commercial products without written approval from Takara Bio Inc. If you require licenses for other use, please contact us by phone at +81 77 565 6973 or from our website at www.takara-bio.com. Your use of this product is also subject to compliance with any applicable licensing requirements described on the product web page. It is your responsibility to review, understand and adhere to any restrictions imposed by such statements. All trademarks are the property of their respective owners. Certain trademarks may not be registered in all jurisdictions.

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