# TaKaRa Taq™ HS Low DNA

**Code No. R090A**  
**Size: 200  $\mu$ l x 5**  
**(for 100 reactions)**

**Shipping at – 20°C**  
**Store at – 20°C**

**Components :**  
**TaKaRa Taq HS Low DNA (2X)**  
**dH<sub>2</sub>O for Low DNA**

**200  $\mu$ l x 5**  
**200  $\mu$ l x 5**

**Lot No.**

**Volume:**

**$\mu$ l**

**Expiration Date:**

## Description :

TaKaRa Taq HS Low DNA is the premix PCR reagent developed by the use of TaKaRa Bio's original technique for purification and DNA inactivation to suppress all detectable DNA in the reagent, derived from host cells and/or contaminated from environment, which can be template on PCR. In addition, this enables you to perform highly sensitive and specific PCR amplification quickly because of the use of modified Taq DNA polymerase having higher extension speed and specificity. This enzyme is designed for hot start PCR using the monoclonal antibody, thus it is possible to prevent non-specific amplification due to mis-priming and/or formation of primer dimer before PCR reaction. Since the antibody is denatured in the initial DNA-denaturation step, the additional activation step is not required. This enzyme has neither 5'-3' exonuclease nor 3'-5' exonuclease activities.

## Origin :

*Escherichia coli* carrying a plasmid that encodes the *Thermus aquaticus* DNA polymerase gene.

## Application :

- For DNA amplification of low copy template by PCR
- For microbial flora analysis

## PCR products :

As most PCR products amplified with TaKaRa Taq HS Low DNA have one A added at 3'-termini, the obtained PCR product can be directly used for cloning into T-vector. Also it is possible to clone the product in blunt-end vectors after blunting and phosphorylation of them.

## PCR test :

Good performance of this product was confirmed by PCR using human genomic DNA (amplified fragment: 2 kb) as templates.

## General reaction mixture for PCR (total 20 $\mu$ l) :

TaKaRa Taq HS Low DNA	10 $\mu$ l
Template	< 100 ng
Primer 1	0.2 $\mu$ M <sup>*1</sup> (final conc.)
Primer 2	0.2 $\mu$ M <sup>*1</sup> (final conc.)
dH <sub>2</sub> O for Low DNA	up to 20 $\mu$ l

Note : Reaction mixture can be set up at room temperature. Be sure to keep all reagents on ice.

\*1 : Adjust the primer concentration to 0.6  $\mu$ M, when you use high speed thermal cycler that ramp temperature speed is over 5°C.

## Primer designing :

Each primers should be designed for T<sub>m</sub> value more than 55°C. The T<sub>m</sub> value should be calculated by below approximation calculation.

$$T_m = 4X \text{ (the number of G or C bases)} + 2X \text{ (the number of A or T bases)} + 35 - 2X \text{ (total number of bases)}$$

Note : In the case of calculating with Nearest Neighbor method, T<sub>m</sub> value of designed primer should be more than 60°C.

## PCR conditions :

2 step PCR

94°C <sup>*2</sup>	5 sec.	} 30 - 35 cycles <sup>*3</sup>
65°C	20 sec./kb	

3 step PCR<sup>\*4</sup>

94°C <sup>*2</sup>	5 sec.	} 30 - 35 cycles <sup>*3</sup>
55°C	1 sec.	
68°C	20 sec./kb	

- \*2 : Be sure to set the denaturation temperature at 94°C. Setting denaturation temperature at over 95°C might cause decrease of reactivity due to inactivation of enzyme.
- \*3 : The setting of cycle number over 40 cycles is not recommended.
- \*4 : In case of using primers which is designed for T<sub>m</sub> value less than 55°C in approximation calculation (or less than 60°C in Nearest Neighbor method), select 3 step PCR reaction.

## NOTICE TO PURCHASER: LIMITED LICENSE

### [P1] PCR Notice

Use of this product is covered by one or more of the following US patents and corresponding patent claims outside the US: 5,789,224, 5,618,711, 6,127,155 and claims outside the US corresponding to expired US Patent No. 5,079,352. The purchase of this product includes a limited, non-transferable immunity from suit under the foregoing patent claims for using only this amount of product for the purchaser's own internal research. No right under any other patent claim, no right to perform any patented method, and no right to perform commercial services of any kind, including without limitation reporting the results of purchaser's activities for a fee or other commercial consideration, is conveyed expressly, by implication, or by estoppel. This product is for research use only. Diagnostic uses under Roche patents require a separate license from Roche. Further information on purchasing licenses may be obtained by contacting the Director of Licensing, Applied Biosystems, 850 Lincoln Centre Drive, Foster City, California 94404, USA.

### [L15] Hot Start PCR

Licensed under U.S. Patent No. 5,338,671 and 5,587,287 and corresponding patents in other countries.

### Note

This product is for research use only. It is not intended for use in therapeutic or diagnostic procedures for humans or animals. Also, do not use this product as food, cosmetic, or household item, etc. Takara products may not be resold or transferred, modified for resale or transfer, or used to manufacture commercial products without written approval from TAKARA BIO INC. If you require licenses for other use, please contact us by phone at +81 77 543 7247 or from our website at [www.takara-bio.com](http://www.takara-bio.com). Your use of this product is also subject to compliance with any applicable licensing requirements described on the product web page. It is your responsibility to review, understand and adhere to any restrictions imposed by such statements. All trademarks are the property of their respective owners. Certain trademarks may not be registered in all jurisdictions.

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