# Yatalase -Plus-

Code No. T030

Size:

2 g

## Description :

This product has complex lytic activities of fungal cell mainly consisting of chitinase, chitinobiase and  $\alpha$ -1,3-glucanase activity.

This product leads to more efficient production of protoplast from filamentous fungi such as *Aspergillus niger* than conventional product.

## Appearance :

Lyophilized powder (containing lactose as the excipient)

## Storage :

4°C, dry condition

## Enzymatic activities :

Chitinase activity :  $\geq 50$  U/g - powder

Chitinobiase activity :  $\geq 500$  U/g - powder

$\alpha$ -1,3-Glucanase activity :  $\geq 20$  U/g - powder

## Unit definition :

[ Chitinase activity ]

1 unit is defined as the amount of enzyme required to release 1  $\mu$  mol of *N*-Acetylglucosamine from 0.5% chitin powder solution in 1 min at 37°C, pH 6.0.

[ Chitinobiase activity ]

1 unit is defined as the amount of enzyme required to release 1  $\mu$  mol of *p*-Nitrophenol from 5 mM *p*-Nitrophenyl-*N*-acetyl- $\beta$ -D-glucosaminide solution in 1 min at 37°C, pH 6.0.

[  $\alpha$ -1,3-Glucanase activity ]

1 unit is defined as the amount of enzyme required to release 1  $\mu$  mol of glucose from  $\alpha$ -1,3-glucan suspension in 1 min at 37°C, pH 6.0.

## Optimum reaction condition :

30°C, pH 6.0

## Quality Control Data :

Please see the Certificate of Analysis (CoA) for each lot. You can download the CoA on Takara Bio website.

## Example of protoplast preparation :

(1) *Aspergillus niger*, *Aspergillus luchuensis*, *Aspergillus sojae*,  
*Aspergillus nidulans*, *Aspergillus kawachii*

|            |                                                                                                                               |
|------------|-------------------------------------------------------------------------------------------------------------------------------|
| Medium     | Dextrin-peptone, 200 mM phosphate Buffer (pH 7.0)                                                                             |
| Culture    | 30°C, 42 hr, Shake culture (Rotary at 140 rpm)                                                                                |
| Conditions | 2% Yatalase -Plus- solution<br>0.8 M NaCl<br>50 mM phosphate Buffer (pH 6.0)<br>30°C Reciprocal shaking (at 60 - 70 rpm) 3 hr |

(2) *Aspergillus oryzae*, *Aspergillus aculeatus*

|            |                 |
|------------|-----------------|
| Medium     | Dextrin-peptone |
| Culture    | Same as in (1)  |
| Conditions | Same as in (1)  |

Manufactured by Ozeki corporation.

## Note

This product is for research use only. It is not intended for use in therapeutic or diagnostic procedures for humans or animals. Also, do not use this product as food, cosmetic, or household item, etc. Takara products may not be resold or transferred, modified for resale or transfer, or used to manufacture commercial products without written approval from Takara Bio Inc.

If you require licenses for other use, please contact us from our website at [www.takarabio.com](http://www.takarabio.com).

Your use of this product is also subject to compliance with any applicable licensing requirements described on the product web page. It is your responsibility to review, understand and adhere to any restrictions imposed by such statements.

All trademarks are the property of their respective owners. Certain trademarks may not be registered in all jurisdictions.

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